Tryptic Soy Agar (Sheep blood 5%)

General purpose medium for the cultivation of a wide variety of microorganisms.

TYPICAL FORMULA

	(g/l)
Tryptone	15.0
Soytone	5.0
Sodium Chloride	5.0
Sheep Blood Defibrinated	50.0 ml
Agar	15.0
Final pH 7.3 ± 0.2	

DESCRIPTION

Tryptic Soy Agar (Sheep blood 5%) is a medium used for the growth of fastidious and non-fastidious organisms. Sheep blood is incorporated to facilitate the growth of some organisms and for the determination of haemolytic reactions.

PRINCIPLE

Tryptone and Sstone provide nitrogen, carbon, sulphur and other essential growth factors. Sodium chloride maintains the osmotic balance of the medium. Sheep blood defibrinated supplies additional growth factors for fastidious microorganisms and allows to evidence the haemolytic reactions. Agar is the solidifying agent.

TECHNIQUE

Inoculate the plates by streaking the sample to examine onto the surface of the medium using a sterile loop in order to isolate single colonies. Plates may be incubated in an atmosphere containing approximately 3-10% carbon dioxide. Incubates at 36 ± 1 °C for 18-24 h.

INTERPRETATION OF RESULTS

Observe for growth and for haemolytic reactions. Four different kinds of haemolysis can be distinguished:

- alpha haemolysis: haemoglobin is reduced to metahaemoglobin in the medium surrounding the colony and this causes a greenish discolouration of the medium;
- beta haemolysis: it is the lysis of erythrocytes which is evident in the bright zone around the colony;
- gamma haemolysis: any destruction of erythrocytes or any change in the medium does not occur;
- alpha-prime haemolysis: a little zone of complete haemolysis surrounded by a partial lysis area is evident.

STORAGE AND TRANSPORT CONDITIONS

2-8°C away from light, until the expiry date on the label. However, our stability studies have shown that the transport at 18-25°C for 4 days, or at 35-39°C for 48 hours, does not alter in any way the performance of the product. Eliminate if signs of deterioration or contamination are evident.

WARNING AND PRECAUTIONS

The product does not contain hazardous substances in concentrations exceeding the limits set by current legislation and therefore is not classified as dangerous. It is nevertheless recommended to consult the safety data sheet for its correct use. The product is designed for *In vitro* diagnostic use and must be used only by properly trained operators.

DISPOSAL OF WASTE

Disposal of waste must be carried out according to the national and local regulations in force.

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Tryptic Soy Agar (Sheep blood 5%)

PRESENTATION

Ready to use plates (90 mm) containing 22 ± 1 ml of medium

STORAGE

2-8°C

PACKAGING

Ref.	Content	Packaging
11037	20 plates	10 plates in thermally soldered film 2 x 10 plates in cardboard box
11037*	100 plates	10 plates in thermally soldered film 10 piles (10 x 10 plates) in cardboard box

pH OF THE MEDIUM

7.3 ± 0.2

USE

Tryptic Soy Agar (Sheep blood 5%) is a general purpose medium used for the cultivation of a wide variety of bacteria and for the determination of haemolytic reactions

TECHNIQUE

Refer to technical sheet of the product

APPEARANCE OF THE MEDIUM

Opaque, cherry red

SHELF LIFE

90 days

QUALITY CONTROL

- Control of general characteristics, label and print
- Sterility control
7 days at $22 \pm 2^\circ\text{C}$, in aerobiosis
7 days at $35 \pm 2^\circ\text{C}$, in aerobiosis
- Microbiological control
Inoculum for productivity: 50-100 CFU
Incubation Conditions: 18-24 hours at $36 \pm 1^\circ\text{C}$

Microorganism		Growth	Haemolysis
<i>Escherichia coli</i>	ATCC® 25922	Good	None
<i>Staphylococcus aureus</i>	ATCC® 25923	Good	Beta
<i>Streptococcus pneumoniae</i>	ATCC® 6305	Good	Alfa
<i>Streptococcus pyogenes</i>	ATCC® 19615	Good	Beta

TABLE OF SYMBOLS

LOT	Batch code	IVD	In vitro Diagnostic Medical Device		Manufacturer		Use by		Fragile, handle with care
REF	Catalogue number		Temperature limitation		Contains sufficient for <n> tests		Caution, consult instructions for use		Do not reuse

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LIOFILCHEM® S.r.l.