STERI TEST VHP MEDIUM (1% Pyr)

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STERI TEST VHP MEDIUM (1% Pyr)

Medium for growth promotion of *Geobacillus stearothermophilus* spores inoculated on biological indicators for monitoring and control of H₂O₂ sterilization processes.

TYPICAL FORMULA (g/L)

Tryptone.....	17.5
Soytone.....	3.0
Sodium Chloride.....	5.0
Glucose	2.5
Dipotassium Phosphate.....	2.5
Bromocresol Purple.....	0.05
Sodium Pyruvate.....	10.0
Final pH	6.6 ± 0.2

DESCRIPTION

STERI TEST VHP MEDIUM (1% Pyr) is a culture medium used for growth promotion of *Geobacillus stearothermophilus* spores inoculated on biological indicators for hydrogen peroxide sterilization processes. It meets the U.S. Pharmacopoeia Revision XXII growth promotion guidelines.

PRINCIPLE

Tryptone and soytone are a source of peptides and aminoacids. Sodium chloride keeps the osmotic balance of the medium. Glucose supplies energy. Dipotassium phosphate is the medium buffer system. Bromocresol purple is the pH indicator. Sodium pyruvate inactivates any hydrogen peroxide residues.

Geobacillus stearothermophilus spores are completely killed off by a sterilization process and therefore, during the following incubation in the **STERI TEST VHP MEDIUM (1% Pyr)** tubes, they are not able to grow and to change the colour of the medium. Otherwise, if the sterilization process is inefficient, the spores partially survive and cause, through glucose fermentation, pH lowering and turn yellow of the medium.

TECHNIQUE

- After submitting the biological indicator to the sterilization process, transfer it aseptically in a **STERI TEST VHP MEDIUM (1% Pyr)** tube.
- Immediately close the tube and incubate it at 55-60°C (131-140°F) for 7 days or for a shorter time validated by user.
- Incubate at the same conditions of time and temperature a biological indicator not submitted to sterilization process in a **STERI TEST VHP MEDIUM (1% Pyr)** tube, belonging to the same batch, as spore growth control (positive control).
- Examine the colour of the medium and interpret results.

INTERPRETATION OF RESULTS

A change of the medium's colour from violet/clear to yellow/turbid indicates spore germination and microbial growth: inefficient sterilization process.

Otherwise, if the colour of the medium remains violet/clear indicates absence of microbial growth: successful sterilization process.

The tube inoculated with the biological indicator not submitted to the sterilization process and used as spore growth control has to turn yellow/turbid after incubation. On the contrary, the test must be repeated after have investigated the cause of the negative result.

STORAGE

2-8°C away from light, until the expiry date on the label or until signs of deterioration or contamination are evident.

WARNING AND PRECAUTIONS

The product is not classified as hazardous by current legislation and does not contain harmful substances in concentration of ≥ 1%. The product must be used only by properly trained operators. Before use ensure the product does not show signs of deterioration and contamination.

DISPOSAL OF WASTE

Disposal of waste must be carried out according to national and local regulations in force.

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STERI TEST VHP MEDIUM (1% Pyr)

PRESENTATION

Ready-to-use glass tubes containing 4 ml of medium.

STORAGE

2-8°C

PACKGING

Code	Content	Packaging
20200	20 tubes x 4 ml	20 tubes in cardboard box

pH OF THE MEDIUM

6.6 ± 0.2

USE

STERI TEST VHP MEDIUM (1% Pyr) is a culture medium used for growth promotion of *Geobacillus stearothermophilus* spores inoculated on biological indicators for hydrogen peroxide sterilization processes. It meets the U.S. Pharmacopoeia Revision XXII growth promotion guidelines.

TECHNIQUE

Refer to technical sheet of the product.

APPEARANCE OF THE MEDIUM

Clear violet medium, without precipitates.

SHELF LIFE

2 years.

QUALITY CONTROL

- Control of general characteristics, label and print.
- Sterility control:
2 days at 20 ± 2°C
1 day at 36 ± 1°C
- Microbiological control:
Inoculum for productivity: 10-100 UFC/ml.
Incubation conditions: 12-18-24 h at 55-60°C.

Microorganism	Growth	Colour
<i>Geobacillus stearothermophilus</i> ATCC 7953	+	yellow

TABLE OF SYMBOLS

LOT	Batch code	 Temperature limitation	 Manufacturer	 Contains sufficient for <n> tests
REF	Catalogue number	 Keep away from heat	 Use by	 Caution, consult accompanying documents

LIOFILCHEM® S.r.l.

Via Scozia, Zona Ind.le - 64026, Roseto degli Abruzzi (TE) - ITALY
Tel +39 0858930745 Fax +39 0858930330 Website: www.liofilchem.net E-mail: liofilchem@liofilchem.net