Acetamide Broth

Liquid medium for detection of *P. aeruginosa* in water samples, according to ISO 16266.

DESCRIPTION

Acetamide Broth is a liquid medium used for confirmation of *Pseudomonas aeruginosa* with the membrane filtration technique.

This medium complies with ISO 16266 for examination of water with low background flora such as waters intended for human consumption or pool waters.

TYPICAL FORMULA*

	(g/l)
Potassium Di-hydrogenphosphate	1.0
Magnesium Sulfate, anhydrous	0.2
Acetamide	2.0
Sodium Chloride	0.2
Sodium Molybdate	0.005
Iron Sulfate, heptahydrate	0.0005
Final pH 7.0 ± 0.5 at 25°C	

*Formula may be adjusted and/or supplemented as required to meet performance specifications;
Grams per litre of purified water.

METHOD PRINCIPLE

Potassium phosphate is the buffering agent. Magnesium sulfate, sodium molybdate and iron sulfate are inorganic salts essential for organisms growth. Acetamide is the sole source of carbon in the medium. Sodium chloride maintains the osmotic balance of the medium. Acetamide is deaminated by gram-negative non-fermenting bacteria to produce ammonia.

PREPARATION

Dehydrated medium Suspend 3.4 g of the powder in 1 liter of distilled or deionized water. Mix well. Heat to boil shaking frequently until completely dissolved. Dispense into tubes or other final containers. Sterilize in autoclave at 121°C for 15 minutes.

TEST PROCEDURE

Filter volumes of diluted or undiluted water sample through 0.45 µm sterile membranes. Place each membrane on a plate of *Pseudomonas* CN agar and incubate at 36 ± 2°C for 40-48 hours. Count all colonies that produce blue/green (pyocyanin) color as confirmed *P. aeruginosa*.

Acetamide Broth is used for confirming fluorescent non-pyocyanin producing colonies and reddish brown pigmented colonies:

- Transfer all colonies to Nutrient agar and incubate at 36 ± 2°C for 20-24 h;
- Inoculate a tube of Acetamide broth with the subculture and incubate at 36 ± 2°C for 20-24 h;
- Add 1-2 drops of Nessler's Reagent and examine for production of ammonia.

INTERPRETING RESULTS

The formation of a yellow to brick red color indicates ammonia production. A negative reaction shows no color change.

Other confirmatory tests, such as Oxidase Test and King's B Medium may be required.

STORAGE

The powder is very hygroscopic, store the powder at 10-30°C, in a dry environment, in its original container tightly closed. Store tubes at 2-8°C away from light. Do not use the product beyond its expiry date on the label or if product shows any evidence of contamination or any sign of deterioration.

SHELF LIFE

Dehydrated medium: 4 years.

Medium in tubes: 1 year.

QUALITY CONTROL

Appearance of dehydrated medium: Free-flowing, homogeneous, white to light beige.

Appearance of prepared medium: Colourless, limpid solution.

Expected cultural response:

Control strain	Inoculum	Incubation	Growth	Acetamide Test
<i>Pseudomonas aeruginosa</i>	≤100 CFU	20-24 h/ 36 ± 2°C	Good	Positive
<i>Pseudomonas aeruginosa</i>				
<i>Pseudomonas aeruginosa</i>				
<i>Escherichia coli</i>				Negative

Please refer to the actual batch related Certificate of Analysis (CoA).

WARNING AND PRECAUTIONS

For professional use only. Operators must be trained and have certain experience. Please read the instructions carefully before using this product. Reliability of assay results cannot be guaranteed if there are any deviations from the instructions in this document.

Consult the Safety Data Sheet (SDS) for information regarding hazards and safe handling practices.

DISPOSAL OF WASTE

Disposal of waste must be carried out according to national and local regulations in force.

BIBLIOGRAPHY

See the references at the end of this document.

TABLE OF SYMBOLS

See the table of symbols at the end of this document.

The product is available in the configurations listed below. There may be additional product ref. numbers as well. For an updated listing of available products, visit [liofilchem.com](https://www.liofilchem.com)

Product	Format	Packaging	Ref.
Acetamide Broth	20x100mm Tube	20 x 5 ml	24154
	16x100mm Plastic Tube	100 x 5 ml	26154P
	Dehydrated medium	500 g	610392

Revision History

Revision	Release Date	Change Summary
1	2024-10-22	Updated layout and content
0.1	2020-05-29	Updated ordering info

This IFU document and the SDS are available from the online Support Center:

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