ESC swab

Demi Fraser Broth

(ISO 18593, ISO 17604)

DESCRIPTION

ESC (Easy Surface Checking) swab - Demi Fraser Broth is a device used for the microbiological monitoring of surfaces by swab sampling method. The medium, also referred to as Half Fraser Broth, is recommended for the detection and enumeration of *Listeria monocytogenes* and *Listeria* spp from food, animal feed, and environmental samples, according to ISO 11290-1 and ISO 11290-2.

CONTENTS OF THE PACKAGES

The product consists of a transparent plastic tube containing liquid medium and a sterile swab attached to the cap.

Each package contains:

- 100 tubes with swab and liquid medium;
- 1 instructions sheet.

METHOD PRINCIPLE

Demi Fraser Broth contains peptic digest of animal tissue, peptic digest of casein, beef and yeast extracts as nitrogen, carbon, vitamin, and mineral sources. Ferric ammonium citrate and esculin are differential agents: all *Listeria* species hydrolyze aesculin to aesculetin which in turn reacts with ferric ions producing a blackening of the medium. Selectivity is provided by the presence of lithium chloride, nalidixic acid, and acriflavine along with an elevated concentration of sodium chloride.

Neutralizing agents, i.e. histidine, lecithin, polysorbate 80 and sodium thiosulfate, are added to the original formula to make this medium suitable for use in areas where residual disinfectants are presumably present.

TEST PROCEDURE

1. Take a tube from the package and unscrew the cap with an attached swab.
2. Press the tip of the swab against the wall of the tube to remove excess liquid.
3. Streak the swab on a surface horizontally and vertically while rotating swab. A Sampling Template 10x10 (ref. 96762) can be used to delimit and sample an area of 100 cm².
4. Return the swab to the tube.
5. Screw the cap and record date and the sampling point.
6. Transport the device to the laboratory and examine according to methods in use.

QUALITY CONTROL

Each batch of ESC swab - Demi Fraser Broth is submitted to quality control using the following strains: *L. monocytogenes* serovar 4b WDCM 00021, *E. coli* WDCM 00013, *E. faecalis* WDCM 00009.

STORAGE

10-25°C away from light, until the expiry date on the label. Eliminate if signs of deterioration or contamination are evident.

WARNING AND PRECAUTIONS

ESC Swab - Demi Fraser Broth does not contain hazardous substances in concentrations exceeding the limits set by current legislation and therefore is not classified as dangerous. It is nevertheless recommended to consult the safety data sheet for its correct use. The product must be used only by properly trained operators.

BIBLIOGRAPHY

- ISO 18593:2018. Microbiology of the food chain - Horizontal method for surface sampling.
- ISO 11290-1:2017. Microbiology of the food chain - Horizontal method for the detection and enumeration of *Listeria monocytogenes* and *Listeria* spp- Part 1: Detection Method.
- ISO 11290-2:2017. Microbiology of the food chain - Horizontal method for the detection and enumeration of *Listeria monocytogenes* and *Listeria* spp- Part 2: Enumeration Method.
- ISO 17604:2015. Microbiology of food the food chain - Carcass sampling for microbiological analysis.
- ISO 11133:2014. Microbiology of food, animal feed and water – Preparation, production, storage and performance testing of culture media.

PRESENTATION	Ref.	Contents
ESC swab - Demi Fraser Broth	85614	100 x 10 ml tubes

TABLE OF SYMBOLS

LOT Batch code	Fragile, handle with care	Manufacturer	Use by
REF Catalogue number	Do not reuse	Contains sufficient for <n> tests	Caution, consult Instruction For Use

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